

HANDLING THE COMFORT CLAMPING DEVICE

Always use both hands to activate the clamping elements simultaneously.

Do not activate both levers (red and green) of one clamping element at the same time!



- ① Lift the red levers for sliding the clamping cover up and down.
Do not touch the green levers when doing that.
- ② When the lid is correctly positioned on the sieve stack, press on the green levers 1 - 2 times to fix the clamping cover on the sieve stack.
Always operate both clamping units simultaneously.
- ③ To loosen the cover after a sieving procedure, lift the red levers and slide the clamping device upwards until the sieve stack can be removed from the sieve shaker. There is no need to take off the device from the support rods completely.

