

# ASSEMBLY OF COMFORT CLAMPING DEVICE



- ① Put both comfort clamping elements on a table, green lever facing down. Place the comfort lid face down on the clamping elements.
- ② Place the O-rings on the cone shaped assembly aid and slide it down into the groove.
- ③ Put the assembly aid in the openings of the lid.
- ④ Place the counterpart of the assembly aid on the cone and slide it down. This presses the O-ring on the clamp and fixes it to the lid. Repeat for the other side.
- ⑤ Slide the clamping cover on the support rods of the sieve shaker, clamping elements facing up.

